07554-80 with pump head 07518-12

OPERATING MANUAL:  
**PUMP DRIVES**

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Edition 02

## SAFETY PRECAUTIONS

**DANGER:** High voltages exist and are accessible in the Console Drive. Use extreme caution when servicing internal components.

### **WARNING: PRODUCT USE LIMITATION**

This product is not designed for, nor intended for use in patient connected applications; including, but not limited to, medical and dental use, and accordingly has not been submitted for FDA approval.

### **Variety of Pump Heads Accepted**

Mount up to 1 (600 rpm) or 2 (200 rpm) MASTERFLEX® L/S® pump heads and all MASTERFLEX-compatible pump heads.

### **Setup and Drive Operation**

1. Mount pump head and load tubing. (See pump head manual.)
2. Adjust flow rate with the 1-turn potentiometer speed control.

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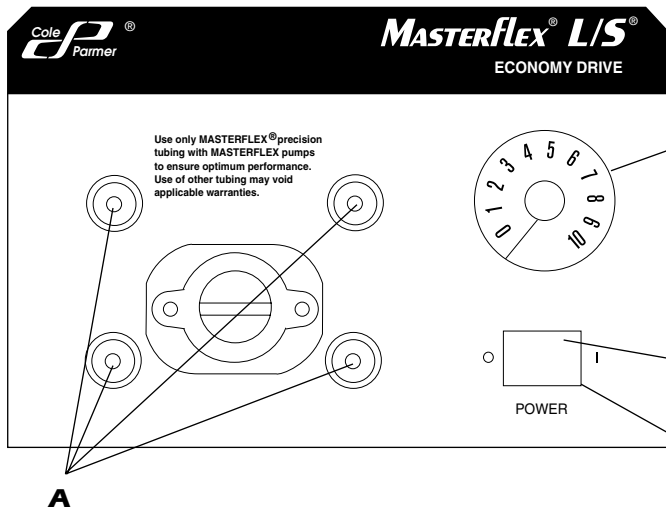
## SAFETY PRECAUTIONS SETUP AND DRIVE OPERATION

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## MOTOR BRUSH CHECK/REPLACEMENT

**NOTE:** Brushes should be checked every 6 months or 2000 operating hours if erratic operation occurs.

**DANGER: High voltages exist and are accessible in the Console Drive. Use extreme caution when servicing internal components.**

1. Place the POWER SWITCH in the off position.
2. Disconnect the AC power input line cord from the AC receptacle.
3. Remove the screws from each side of the housing and lift off the housing.
4. Carefully unscrew each brush cap. Withdraw the brush, and examine it for wear.

**NOTE:** Replace both brushes, if either brush is less than 7.6 mm (0.300 in) long from base to point.

5. Insert brushes and install brush cap.
6. Install housing and secure with the screws on each side.
7. Connect the AC power input line cord to the AC receptacle.

## Fuses

**NOTE:** These drives have current limiting which cause drive to stop under overload conditions. If drive does not start return to dealer for servicing.

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## MOTOR BRUSH CHECK/REPLACEMENT CLEANING

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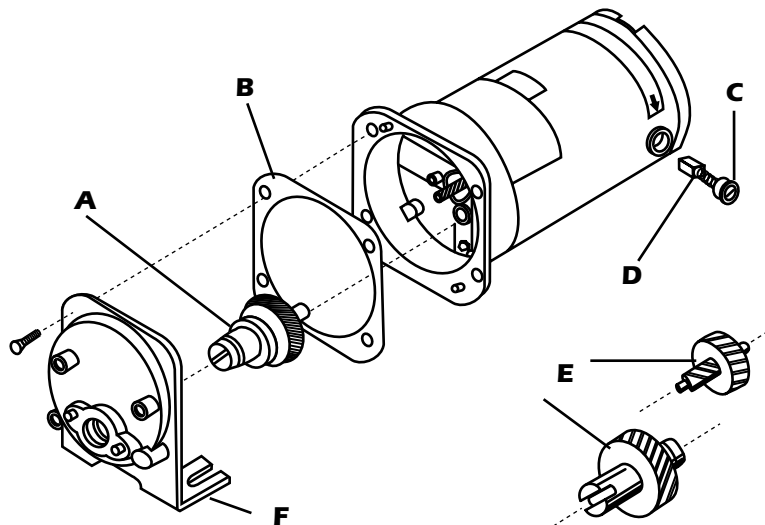
<b>Power Output:</b>	<b>37 W (0.05 hp)</b>
<b>Maximum Current:</b>	
115 V Units:	1.5A, shorted output conditions
230 V Units:	0.9A, shorted output conditions
<b>Speed Regulation:</b>	
Line:	±2%
Load:	±3%
Temperature Drift:	±10%
<b>Enclosure Rating:</b>	<b>IP22 per IEC529</b>
<b>Humidity (non-condensing):</b>	<b>10% to 90%</b>
<b>Operating Temperature:</b>	<b>0° to 40° C</b>
<b>Storage Temperature:</b>	<b>-45° to 60° C</b>
<b>Chemical Resistance:</b>	Exposed material is painted steel housing
<b>Altitude:</b>	Less than 2000 m
<b>Display:</b>	Green LED
<b>Dimensions (L × W × H):</b>	220 mm × 180 mm × 135 mm (8 <sup>11</sup> / <sub>16</sub> in × 7 <sup>1</sup> / <sub>8</sub> in × 5 <sup>5</sup> / <sub>16</sub> in)
<b>Weight:</b>	4.1 kg (9 pounds)
<b>Compliance:</b>	115V: UL508, CSA C22.2, No. 14-M91 230V (For CE Mark): EN61010-1/A2: 1995 (EU Low Voltage Directive) and EN61326-1/A1: 1998 (EU EMC Directive)
<b>Pollution Degree:</b>	Pollution Degree 2 per IEC 664 (Indoor usage—lab, office)
<b>Installation Category:</b>	Installation Category II per IEC 664 (Local level—appliances, Portable Equipment, etc.)

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## EU Declaration of Conformity

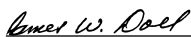
**Name of Apparatus:** MASTERFLEX® Economy Analog Console L/S® Pump Drives  
**Model Numbers:** 7554-85 & -95, 77910-25  
**Description of Apparatus:** Variable speed, peristaltic pump motor drive. Used with pump head and peristaltic tubing to pump fluids.  
 7554-85 (200 rpm) & -95 (600 rpm);  
 77910-25; (Bundled 7554-95)

Barnant Company declares that the above models are in conformity to the following harmonized standards and directives

Applicable Directives	Applicable Specifications	Manufacturer's Report Number
73/23/EEC 93/68/EEC	EN61010-1/A2:1995	TR9651
89/336/EEC 92/31/EEC 93/68/EEC	EN61326-1/A1:1998	TR9652

The last two digits of the year in which the current configuration of the above models were assessed per the Low Voltage Directive is :00.

**Manufacturer:** Barnant Company Division  
 Cole-Parmer Instrument Company  
 28W092 Commercial Avenue  
 Barrington, IL 60010-2392  
 USA  
 Tel: 847-381-7050

**Manufacturer's Signature:**  17 July, 2000  
 James W. Doll Date  
 Vice President, Engineering

## WARRANTY

Use only MASTERFLEX precision tubing with MASTERFLEX pumps to ensure optimum performance. Use of other tubing may void applicable warranties.

The Manufacturer warrants this product to be free from significant deviations from published specifications. If repair or adjustment is necessary within the warranty period, the problem will be corrected at no charge if it is not due to misuse or abuse on your part, as determined by the Manufacturer. Repair costs outside the warranty period, or those resulting from product misuse or abuse, may be invoiced to you.

The warranty period for this product is noted on the Warranty Card.

## PRODUCT RETURN

To limit charges and delays, contact the seller or Manufacturer for authorization and shipping instructions before returning the product, either within or outside of the warranty period. When returning the product, please state the reason for the return. For your protection, pack the product carefully and insure it against possible damage or loss. Any damages resulting from improper packaging are your responsibility.

## TECHNICAL ASSISTANCE

If you have any questions about the use of this product, contact the Manufacturer or authorized seller.

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